

# Surface RNase Decontamination



## DECON-R™

Used to remove traces of RNASE and virus contamination from laboratory platforms, tiles, work bench surfaces, laboratory apparatus, equipments and glass/plastic surfaces

### Instructions for Use

**On Work Surfaces:** Apply **DECON-R™** solution directly on the surfaces to be cleaned, wipe the work surface thoroughly using a clean paper/cloth towel, rinse with water and wipe dry using another clean paper/cloth towel.

**For Laboratory Apparatus:** Apply liberally **DECON-R™** solution to a clean paper/cloth towel, wipe the surface of the apparatus with this towel. Then rinse with water and wipe dry using another clean paper/cloth towel. Alternatively, small parts of the apparatus may be cleaned by soaking these parts in **DECON-R™** solution for minimum 10 minutes, followed by rinsing with water and then wipe dry using another clean paper/cloth towel.

**For Plastic/Glass Equipment Surfaces/Centrifuges:** Rub enough **DECON-R™** solution on to the equipment surfaces, such that the entire surface gets coated with the solution. Rinse it thoroughly with distilled water at least two times, then wipe dry it completely. It is important to wear appropriate protective equipments such as safety glass, gloves or protective clothing while handling chemical.

### Presentation

REF	1108092000
Σ	10 x 200 ml